

Student Slide Coverslips MS-SLIDCV

Cleaning Instructions

These value priced student grade coverslips are not pre-cleaned. Although not usually necessary, cleaning with a light soap and water may be required for best optical quality. Follow these steps if cleaning is required:

1. Add 1-2 drops dishwashing liquid detergent to a bowl of lukewarm water. Swish the soap around in the bowl to produce suds.
2. Add a few slide coverslips to the dish and very gently rub each coverslip between your thumb and forefinger.
3. Remove the coverslips from the bowl and rinse in a bowl of cold water.
4. Carefully pat dry each coverslip with a soft cloth or paper towel to avoid water spots.

CAUTION: Slide coverslips are very thin and can break easily. Use care when following this procedure to avoid breaking the coverslips and cutting yourself on the broken glass.